

automatic serial determination of up to 100 samples rapid-clean-precise



N-REALYZER

Elemental analysis for the determination of nitrogen

Technical data N-Realyzer

Data	N-Realyzer
Sample size	up to 1000 mg
Sample capacity	· •
	3 - 5 min, depending on the sample type and sample size
Recovery rate	
Detection limits with Helium	0,003 - 50 mg N absolut
Standard deviation	< 0,5 % (with standard sample preparation)
Operation	via computer with REAL-OS
Temperature range Combustion furnace	400 - 1.100 °C
Temperature range Reduction furnace	400 - 1.100 °C
Temperature range Desorption furnace	50 - 400 °C
Required gases and quality	Helium, quality grade 5.0 (99,999 %)
	Oxygen, quality grade 5,0 (99.999 %)
	Compressed air (100 % oil- and water free)
Gas installation	ressure regulators for high purity gas; 1/4" cylindric thread (female)
Inlet pressure helium	3 - 6 bar
Inlet pressure oxygen	3 - 6 bar
Inlet pressure compr. air / nitrogen	7 - 8 bar
Ambient temperature	15 °C < t < 35 °C
Nominal voltage	100 - 240 VAC
Frequency	50/60 Hz
Dimensions Device closed W / D / H	
Dimensions Device opened* W / D / H	600 x 550 x 1200 mm
3	ca. 130 kg
Nominal wattage	
Current	max. 14 A (6 A bei 230 V, 14 A bei 100 V)
Wanted digital balance	Precision min. 0,1 mg

^{*} Upper protective cover open



ORDER INFORMATION

When ordering the apparatus, please note that you must order the correspondin REAL-OS control unit for the intended destination country in addition to the 14-4000 base unit. The base unit can be operated for all voltages from 100 - 240 volts. The respective control units are configured for the corresponding target markets with suitable connection cables.

N-Realyzer Base unit



Order No.	Туре	Description
14-4000	REAL-N	N-Realyzer basic unit, complete with: - Autosampler for the automatic supply of up to 100 samples - Starter kit incl. all consumables for 1,000 samples [absorber, CO ₂ adsorber, copper, ash insert, etc.] - Accessories and tools for sample handling - Connection lines for gas and compressed air installation

REAL-OS Control Unit 230 V (EUR/INT)






Plug type F 230 V

Order No.	Туре	Description
14-4002	CUN-02	REAL-OS Control Unit for N-Realyzer, for target markets Europe and International with plug connection type Typ F, 230 V - complete with: - 14-4545 Mini PC with REAL-OS control software and power supply unit - 14-4549 Ethernet cable, 2 m (2 pieces) - 10-0057 Mains cable AKE 3x1,0 mm², 1,5 m Note: Monitor, PC keyboard and mouse are not included in the scope of delivery and must be provided on site*

REAL-OS Control Unit 220 V (US/CAN/CSA)

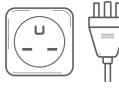


Plug type B, NEMA 6-15 220 V

Order No.	Туре	Description
14-4003	CUN-03	REAL-OS Control Unit for N-Realyzer, for target markets USA, Canada, South America with plug connection type NEMA 6-15, 230 V - complete with: - 14-4545 Mini PC with REAL-OS control software and power supply unit - 14-4549 Ethernet cable, 2 m (2 pieces) - 1004799 US device connection cable NEMA 6-15 with plug, 2.5 m - 14-4559 Travel plug N-Realyzer Control Unit Note: Monitor, PC keyboard and mouse are not included in the scope of delivery and must be provided on site*



REAL-OS Control Unit 115 V (US/JPN/CSA)

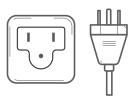

Plug type B, NEMA 5-15 115 V Attention, only permitted with protective conductor!

Onder Ne	T	December 1
Order No.	Туре	Description
14-4004	CUN-04	REAL-OS Control Unit for N-Realyzer, for target markets USA, Japan, South America with plug connection type B, NEMA 5-15, 115 V - complete with: - 14-4545 Mini PC with REAL-OS control software and power supply unit - 14-4549 Ethernet cable, 2 m (2 pieces) - 1007344 15A mains cable, with NEMA 5-15, 115V [with IEC plug] - 14-4559 Travel plug N-Realyzer Control Unit Note: Monitor, PC keyboard and mouse are not included in the scope of delivery and must be provided on site*

REAL-OS Control Unit 240 V (GB)





Plug type G 240 V

Order No.	Туре	Description
14-4005	CUN-05	REAL-OS Control Unit for N-Realyzer, for target markets Great Britain with plug connection Type G 1-13 A, 240 V - complete with: - 14-4545 Mini PC with REAL-OS control software and power supply unit - 14-4549 Ethernet cable, 2 m (2 pieces) - 10-0050 Mains cable, 1.5 m with English plug plug Note: Monitor, PC keyboard and mouse are not included in the scope of delivery and must be provided on site*

^{*}Minimum monitor requirements: Full HD resolution 1080 p (1920 x 1080), 24" screen diagonal, recommended: HDMI interface, Vesa mount compatible, 75 Hz

^{*}Minimum keyboard and mouse requirements: 100% keyboard with Numpad, wireless



ACCESSORIES AND CONSUMABLES

The best analytical results can only be achieved with original accessories from C. Gerhardt. These have been specially developed and tested for use with the N-Realyzer in our in-house application laboratory.

Set of Consumables - DumaSet (with high-performance reactor)



Order No.	Туре	Description
14-0194	DumaSet	N-Realyzer - Compilation of important consumables, consisting of: 1 x 14-0244 DumaReact, combustion reactor, filled with HT and LT 2 x DumaTube, quartz tube for reactor 1 x DumaCollect, ash insert (unit with 10 pcs.)
		10 x DumaFoil, Tin foil, conditioned, (10 boxes á 100 pcs.)

Reactor - DumaReact



Order No.	Туре	Description
14-0244	DumaReact	Prepacked combustion reactor,
		filled with HT- and LT-catalyst, 1 pc.

Quartz tube - DumaTube



Order No.	Туре	Description
14-0203	DumaTube	Quartz tube for reactor, OD 26mm, L 450mm, 2mm wall thickness, 1 pc

Copper - DumaCop



Order No.	Туре	Description
14-0007	DumaCop	Copper for reduction, 10 bottles á 225 g (two bottles are enough for one filling for quartz tube no. 14-0203)
14-0002	DumaCop	Copper for reduction, 20 bottles á 225 g (two botles are enough for one filling for quartz tube no. 14-0203)

HT wool - DumaPad (for quartz tube and combustion reactor)



Order No.	Туре	Description
14-0225	DumaPad	HT Wool pads combibox, content: 30 pcs. small + 30 pcs. large
14-0274	DumaPad	HT Wool pads small , content: 60 pcs.
14-0275	DumaPad	HT Wool pads large , content: 60 pcs.



Special tool for N-Realyzer



Order No.	Туре	Description
14-0064		Tool for taking out reactor tubes,1 pc.

Removing tool



Order No.	Туре	Description
14-4511		Removing tool for ash insert

Extractor



Order No.	Туре	Description
14-0068		Extractor for ash insert, 1 pc. For removing and inserting hot ash inserts

Tin foil - DumaFoil



Order No.	Туре	Description
14-0017	DumaFoil	DumaFoil, tin foil 3.5 x 3.5cm, 100 pcs.
14-0417	DumaFoil	DumaFoil XL, tin foil 4.3 x 4.3 cm, specially designed for weighing in larger samples, 100 pcs.

Ash insert - DumaCollect



Order No.	Туре	Description
14-0015	DumaCollect	Ash insert, ceramics, with bottom, for single use
		L 100 mm, package with 10 pcs.



Condensate collection glass



Order No.	Туре	Description
14-4331		Condensate collection glass, square 50 ml

Standard - DumaEDTA



Order No.	Туре	Description
14-0032	DumaEDTA	EDTA Calibration Standard, 10g

Absorbent - DumaSorb



Order No.	Туре	Description
14-0022	DumaSorb	Absorbent for liquid samples, content 25 g
14-0295	Super- absorber	Absorbent, special, for low salt- and low fat liquid samples of up to max. 400 mg, content 50 g

Adsorber - DumaCO₂



Order No.	Туре	Description
14-4085	DumaCO ₂	REAL-N Spiral adsorber for CO ₂ , completely filled with adsorbent, gas-tight seal incl. GL 14 screw caps, package with 7 pieces
1003687		REAL-N Screw cap GL 14, closed

Probentransfereinheit



Order No.	Туре	Description
14-4529		REAL-N Sample transfer unit for collecting and safely transferring up to 100 samples into N-Realyzer analysis
14 4527		,



Encapsulating device



Order No.	Туре	Description
14-0025		Encapsulating device for samples, 3-part, for sample magazine 64-place

Accessories for water adsorption trap

Order No.	Туре	Description
14-0219		Adsorbent for water adsorption trap 14-0061, Content 100 g, 1 bottle
14-0217		Glass part L = 150 mm for water absorption trap 14-0061
14-0243		Covering wool for water absorption trap, 5g
14-0081		Silicone sealing for water absorption trap, 2 pcs









Rack versions

Order No.	Туре	Description
14-0672		Reactor rack - for the safe storage of up to 4 reactors (DumaReact)
14-0076		Metal stand for supporting reactor tubes







Filling funnel



Order No.	Туре	Description
14-0218		Funnel for the filling of the reactors and traps, 1 pc.

Sampling devices for solid & liquid samples

Order No.	Туре	Description
14-0073		Spatula for solid samples, 1 pc.
14-0095		Syringe, 1 ml for liquid samples, 1 pc.





Rod for filling



Order No.	Туре	Description
14-0065		Rod for reactor tube filling 1 pc

Tweezer



Order No.	Туре	Description
14-0072		Tweezer with curved tip, 1 pc.

End caps for reactor



Order No.	Туре	Description
14-0273		End caps for reactor small/large, content: - 10 x end caps small - 10 x end caps big



Reactor connection, complete



Order No.	Туре	Description
14-4518		Reactor connection, completely assembled with: Teflon tube 14-4436, glass connector 14-4435 and o-ring seal 14-4437

Spare parts for reactor connection

Order No.	Туре	Description
14-4436		Teflon tube for reactor connection, inner Ø 4 mm, complete with nuts and ferrules
14-4435		Glass connector for reactor connection
14-4437		O-ring seal green FFPM, green 18 x 4 mm, pack of 4 pcs.







Leakage detector



Order No.	Туре	Description
14-0008	LD239	Leak detector for detecting leaks in the apparatus, complete with: - USB charging cable

Sealings for quartz tubes

Order No.	Туре	Description
14-0020		Sealings for quartz tube top, black, Ø 26 mm, 4 pcs.
14-4437		Sealings for quartz tube bottom, green, Ø 18 mm, 4 pcs.







Adapter for gas pipes

Order No.	Туре	Description
14-0126		Adapter 1/4", for PU gas pipe 4 mm
14-0125		Adapter 1/4", for gas pipe 2 mm
14-0128		Adapter 3/8", for PU gas pipe 4 mm
14-0127		Adapter 3/8", for gas pipe 2 mm






Spare parts for gas pipes

Order No.	Туре	Description
14-0069		Sealing cone for 4 x 0.75 mm for PU tube, 6 pcs.
14-0070		Nut M6 x 0.75 mm, for 2 mm stainless steel pipe, 10 pcs.
14-0071		Sealing cone for 2 mm stainless steel pipe, 10 pcs.





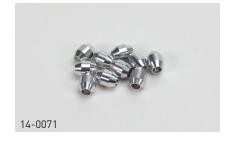



SUPPORT AND SERVICE

In addition to its high quality instruments, C. Gerhardt offers comprehensive services to maintain the functionality and effeciency of your equipment.

INSTALLATION / EQUIPMENT QUALIFICATION

Selection	Туре	Description
99-0204	ABI-DT	Installation and Instruction N-Realyzer - Installation and connection of the unit - functional check - operating instructions
99-0212	GQF-DT	Equipment Qualification (IQ/OQ) N-Realyzer - Installation of the unit and instruction - Installation Qualification IQ and Operational Qualification OQ, Guidance for Performance Qualification PQ - Training of laboratory personnel with certificate - Individual training certificates, if reqested

SERVICING AND MAINTENANCE CONCEPTS

For detailed information on servicing and maintenance concepts, please contact your local authorised C. Gerhardt representative.

WORKSHOPS / SEMINARS / APPLICATIONS TRAINING

For detailed information on trainings and workshops, please contact your local authorised C. Gerhardt representative.

Product data sheet N-Realyzer



TABLE OF CONFIGURATION

Please arrange the required system

N-F			

Quantity	Order No.	Туре	Description

CONSUMABLES / ACCESSORIES

Quantity	Order No.	Туре	Description

SUPPORT AND SERVICE

Quantity	Order No.	Туре	Description

Send offer to:

Customer No.	
Company	
Name	
Address	
Tel.	
E-mail	
Web	



C. Gerhardt GmbH & Co. KG Cäsariusstraße 97 · 53639 Königswinter, Germany Tel.: +49 (0) 2223 2999-0 · www.gerhardt.de

